













Heat Snapshots

Which place has the most heat?

<p>Hold the probe on a radiator</p>  <p>winter only!</p>	<p>Count to 10 then press Green</p> 
<p>Hold the probe in cold water</p>  <p>C</p>	<p>Count to 10 then press Green</p> 
<p>Hold the probe in hot water</p>  <p>H</p>	<p>Count to 10 then press Green</p> 
<p>Hold the probe in a fridge</p> 	<p>Count to 10 then press Green</p> 
<p>Hold the probe in a hand</p> 	<p>Count to 10 then press Green</p> 
<p>Your idea</p>	<p>Count to 10 then press Green</p> 
<p>Your idea</p>	<p>Count to 10 then press Green</p> 

Teacher Notes:

Do not plug the datalogger into the computer.

- Switch on the datalogger with the Green start button.
- Press the blue menu button twice until it says "Press Green to start snapshot >"
- Press the Green button. You should see it say "Logging started" and then the display returns to normal.
- Press Green button each time you take a snapshot. Press the Red button when you have finished – it should say "Logging stopped".
- Plug the datalogger into the computer, open the LogIT software and press Fetch to see the bar graph.