












# Light Snapshots

Which place has the most light?

Point the sensor to a window 	Press Green 
Point the sensor to a computer monitor 	Press Green 
Point the sensor to a torch 	Press Green 
Point the sensor under a table 	Press Green 
Point the sensor outside 	Press Green 
Your idea	Press Green 
Your idea	Press Green 

Teacher Notes:

*Do not plug the datalogger into the computer.*

- Switch on the datalogger with the Green start button.
- Press the blue menu button twice until it says "Press Green to start snapshot >"
- Press the Green button. You should see it say "Logging started" and then the display returns to normal
- Press Green button each time you take a snapshot. Press the Red button when you have finished – it should say "Logging stopped".
- Plug the datalogger into the computer, open the LogIT software and press Fetch to see the bar graph.